




Shedding New Light On MICROSCOPY

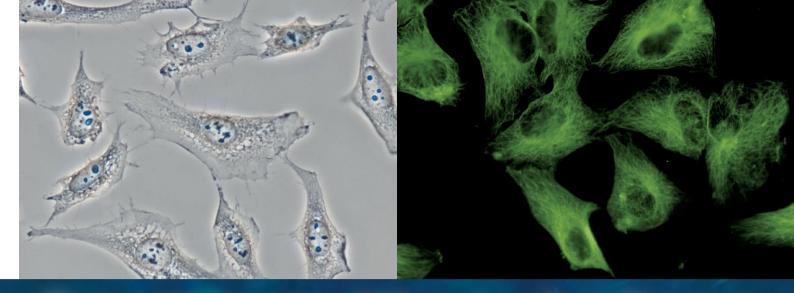

Do more than before

Easy to work with

Simple operations

Control buttons on the Ts2R microscope are intuitively located for a streamlined workflow. The on/off and diascopic/epi-fluorescence

switching controls are located on the front panel while buttons pertaining to either diascopic or epi-fluorescence control are zoned to the left and right sides of the microscope body, respectively.



Mechanical stage

The Ts2R can be configured with the high-grade rectangular mechanical stage. This stage provides a long travel stroke, enabling users to observe an entire well plate from end-to-end. You can also set limits to the travel stroke (three-way) to match frequently used samples or vessels. The stage handle is offered in two lengths, long or short, to further accommodate the user's needs.



Ergonomic stage design for improved workflow

The Ts2R's stage height has been lowered by approximately 30% compared to the standard research microscope*, thereby ensuring a comfortable hand position during repetitive operation and sample exchange. Moreover by lowering the support columns and positioning the camera port on the side of the tube, sample visibility is improved.

*Comparison with Nikon's Inverted Research Microscope ECLIPSE Ti2.



Versatile applications with LED illuminators

The high-intensity LED light source enables you to perform a wide variety of observation methods similar to the full-size, inverted research microscope ECLIPSE Ti2. The Ts2R-FL model, which offers epi-fluorescence observation, provides four fluorescence channels and 8 different wavelengths to choose from.

	Ts2R	Ts2R-FL
DIA	Brightfield	Emboss Contrast Phase Contrast NAMC (Nikon Advanced Modulation Contrast)
FL	_	Epi-fluorescence





A compact inverted microscope for your basic research needs

Compact body

Compact body for streamlined workflow

Illumination modules including the epi-fluorescence light source have been seamlessly incorporated into the microscope main body, resulting in a compact and simple design form that's also durable. The compact structure is also vibration-resistant to provide highly stable sample observations.



Easily fits inside laminar flow hoods

The low stage and side-port camera position reduce user fatigue from repetitive stage manipulation and provides clear visibility of the stage

and sample even with the hood sash lowered.
Additionally, by rotating the eyepiece tube 180° and fastening it in position, it is possible to have the microscope completely within the hood.



High precision and quality

Advanced-optical performance

The Ts2R is compatible with Nikon's acclaimed CFI60 objective lenses which provide high numerical apertures and long working distances to deliver stunningly clear images.

High performance and quality optical accessories

Optical accessories achieve the same performance level as Nikon's inverted research microscope ECLIPSE Ti, providing exceptionally clear, sharp images.

Emboss contrast image (Top left): Two-cell stages derived from mesophyll protoplasts from *Nicotiana benthamiana* after 7 days of culture in liquid medium. (Objective: CFI S Plan Fluor ELWD 20XC)

Photo courtesy of: Dr. Jutta Schulze, Institute of Plant Biology, Braunschweig University of Technology





Do more than before — DIA




Diascopic observations with enhanced quality


High-intensity LED Eco-illumination

Nikon's LED Eco-illumination is environmentally friendly with its low power consumption and yet provides extremely bright illumination, suitable for phase contrast and DIC imaging. The built-in fly-eye lens ensures uniform brightness across the entire field of view. Furthermore, LED excitation has no unwanted UV component, thereby eliminating UV-mediated cell damage and improving cell survival rates during long-term imaging.

Phase contrast observation

Phase contrast is an optical contrasting technique that typically utilizes a phase contrast objective lens and condenser annulus. The use of a high-intensity LED light source results in clear images even at high magnifications.

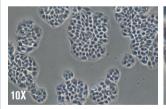

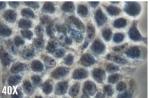

Apodized Phase Contrast (APC) observation

APC observation is a type of phase contrast microscopy which minimizes unwanted halos in thick specimens. For example, APC technique provides clearer details of thick samples such as dividing cells.

Nikon Advanced Modulation Contrast (NAMC)

NAMC provides high relief, DIC-like images of samples on plated on plastic dishes, which is not possible with DIC observation. Ts2R provides high-quality NAMC images like Nikon's inverted research microscope, ECLIPSE Ti2.



Application



New contrasting technique, "Emboss Contrast"

Nikon's new contrasting technique is compatible with both plastic and glass culture dishes. Unlike phase contrast or NAMC, Emboss Contrast does not require special objective lenses and therefore has minimal effect on epi-fluorescence observation. Emboss Contrast allows thick samples such as embryos to be easily observed in pseudo-three-dimensional image with great clarity.

Image courtesy of Hideaki Watanabe, Ph.D. and Hisataka Hasegawa, Ph.D.

Spindle observation


Accurate observation of spindle bodies is easily attained with the Ts2R. The system offers finely detailed work without damaging the spindle body.



Differential Interference Contrast (DIC) observation

DIC provides high-resolution pseudo-three dimensional images that have a shadow-cast appearance. New high-intensity LED illumination results in vivid DIC images even at high magnifications.



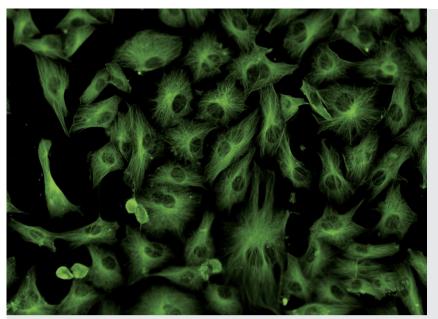




Do more than before — FL



Fluorescence images with uniform bright illumination


Application

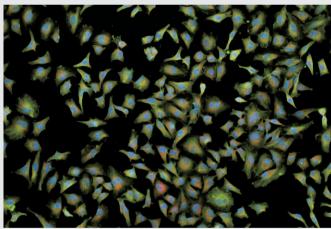

Overlapping image with three colors with use of Imaging Software NIS-Elements

Multicolor fluorescence observation

Using four different LEDs, multicolor fluorescence observation can be easily and efficiently achieved.

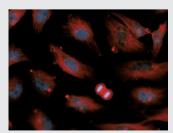

High signal-to-noise fluorescence imaging

Noise Terminator helps to capture vivid images.



Clear, vivid observation across the entire field of view

The fly-eye lens delivers uniform brightness to the entire field of view.

Accurately reproduce illumination power every time

The illumination power previously defined by the user is replicated when the same wavelength is used again, thus eliminating the need


for manual adjustment of light intensity when switching between wavelengths.



High S/N epi-florescence observation in bright rooms

The new Contrast Shield accessory (optional) blocks room light, providing an easy and cost-effective method for achieving high signal-to-noise fluorescence observation in a brightly lit laboratory.



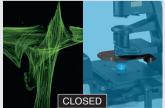

Fluorescent LED light source

The D-LEDI Fluorescence LED Illumination system can be attached for simultaneous fluorescent and phase contrast observation or fluorescent and differential interference contrast observation.



Accessories

ThermoPlate® TPi-TCSX

ThermoPlate® TPi-TCSX provides accurate and stable temperature control for the specimen from room temperature to 50 degrees Celsius. Proprietary treatment methods ensure that the glass surface of the Termo

Manufacturer: TOKAI HIT Co., Ltd.

Plate is breakage-free.







Stage-top incubator -----

A stage-top incubation chamber can be utilized to accurately control temperature, humidity and CO2 levels to maintain optimal cell health during long-term observation.

Manufacturer: TOKAI HIT Co.,Ltd.



This compact manipulation system features a suspension-type, soft-touch joy-stick. The hydraulic remote controls enable smooth, movement-free manipulation, minimizing needle deflection. Users can seamlessly switch between coarse and fine motion. Additionally, indicators on the coarse control mechanisms aid needle adjustments.

Manufacturer: NARISHIGE LIFEMED CO., LTD.



Cameras for microscopes

All cameras of the digital sight series can be directly connected to a PC via a fast USB3.0 interface.

Monochrome Microscope camera

16.25 megapixel

Monochrome

Cooled

Capable of expressing images as

is, this microscope digital camera

offers high-sensitivity, and low noise

DS-Qi2

*The optional camera port is required to attach the digital camera to the microscope.

F-mount CMOS Camera

Microscope camera DS-Ri2



Max recordable pixels



Capable of expressing images as is, this microscope digital camera

offers high resolution, superior color

reproduction and fast frame rate. imaging. Frame rate

6 fps (4908 \times 3264), 45 fps (1636 \times 1088) 4908 × 3264

C-mount CMOS Camera Microscope camera DS-Fi3





A high-definition 5.9-megapixel color CMOS image sensor captures fine-textured images in faithful color. For image acquisition, NIS-Elements imaging software is required.

15 fps (2880 \times 2048), 30 fps (1440 \times 1024) 2880×2048





Simply installing NIS-Elements L on a tablet PC enables setting and control of DS-Fi3/DS-Ri2 microscope cameras, live image display, and image acquisition.

*For information about compatible tablet PCs, contact Nikon.



<u>Syst</u>em diagram

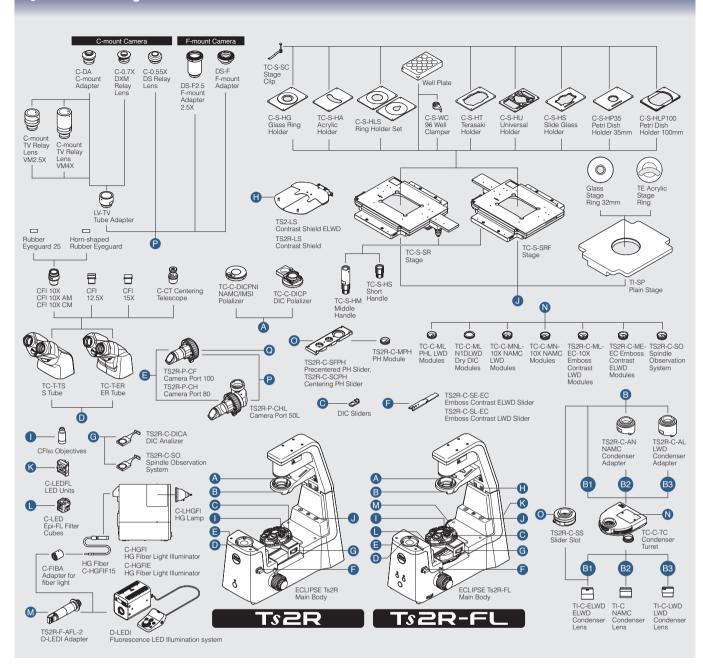

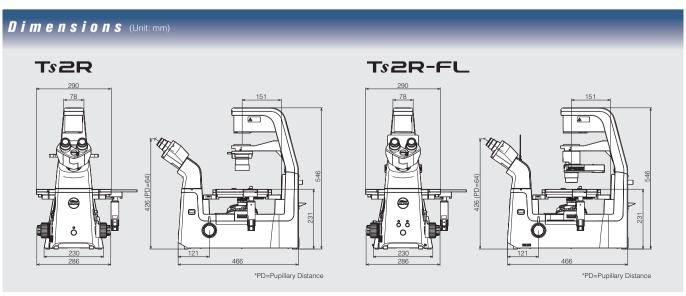

Ts2R / Ts2R-FL Specifications				
	T ₅ 2R	Ts2R-FL		
Optical System	CFIso Infinity Optical System			
Observation method	Brightfield, Apodized Phase Contrast*1, Phase Contrast, Nikon Advanced Modulation Contrast*2, DIC, Emboss Contrast*3, Spindle Observation	Brightfield, Apodized Phase Contrast* ¹ , Phase Contrast, Nikon Advanced Modulation Contrast* ² , DIC, Emboss Contrast* ³ , Epi-Fluorescence, Spindle Observation		
Illumination Diascopic illumination	High luminescent white LED illuminator (Eco-illumination), Built-in Fly eye lens			
Episcopic illumination	_	LED illuminator, built-in Fly eye lens, Can be configured with up to 4 different fluorescence LED units; available wavelengths: 385, 455, 470, 505, 525, 560, 590, 625 nm		
Tube	Binocular tube: Inclination: 35 degree Frgonomic tube: Inclination: 15-45 degree, Siedentopf type, Pupillary distance: 50-75 mm, Attachable camera port			
Eyepiece(F.O.V.)	10X (22), 12.5X (16), 15X (14.5)			
Camera port (Eyepiece: Port)	• TS2R-P-CF 100:0 / 0:100 • TS2R-P-CH 100:0 / 20:80 • TS2R-P-CHL 100:0 / 50:50			
Focusing	Via nosepiece up/down movement, Stroke (manual): Up 8 mm, down 3 mm Coarse stroke: 5.0 mm per rotation, Fine stroke: 0.1 mm per rotation, Coarse motion torque adjustable, Refocusing mechanism mounted			
Nosepiece	Sextuple nosepiece, With DIC prism slots			
Condenser	Condenser turret, mount up to 7 modules: Phase Contrast, DIC, NAMC, IMSI, Emboss Contrast and ND for Bright Field Use with any one of ELWD condenser lens, LWD condenser lens and NAMC condenser lens			
Slider	Precentered or Centering PH Slider, 10X, 20X, 40X Objectives available for phase contrast Emboss Contrast sliders (eyepiece-tube-side slider must be mounted), 10X, 20X, 40X, 60X objectives available for Emboss Contrast			
Stage	Plain Stage, Stage Size 260(X)×300(Y) mm with 2 types of Stage Ring Rectangular Mechanical Stage Stroke: 114(X)×73(Y) mm, Adjustable XY stroke limit, Accepts 8 types of micro-testplate, well clamper and stage clip			
Holder	C-S-HP35 Petridish Holder 35 mm C-S-HLP100 Petridish Holder 100 mm C-S-HT Terasaki Holder for Terasaki holder and Ø65 dish C-S-HS Slide Glass Holder for glass slides, Ø54 dish and hemocytometer C-S-HU Universal Holder for Terasaki plate holder, glass slide, Ø35-65 dish and hemocytometer C-S-HG Glass Ring Holder C-S-HLS Ring Holder Set TC-S-HA Acrylic Holder			
Epi Fluorescence attachment	_	Epi-fluorescence filter turret (with main body), Filter cubes with noise terminator mechanism Configure with up to 4 Epi-fluorescence filter cubes, one position is used during bright-field observation, Attachable Contrast Shield (optional; LWD,ELWD) In combination with the dedicated adapter, D-LEDI Fluorescence LED Illumination system can be used.		
Dimensions	286(W)×466(D)×542(H) mm	286(W)×466(D)×542(H) mm		
Weight (approx.)	17 kg	18 kg		
Rated Voltage/Electric Current	100V-240 V, Less than 0.65 A			
Power Consumption	30W			

^{*1} APC (Apodized Phase Contrast) is a type of phase contrast observation with reduced halo, thanks to Nikon's unique lens coating.

Related Products

ECLIPSE TS2/TS2-FL

Fits in Every Laboratory - Simple to Use & Compact

The new Inverted Routine Microscope ECLIPSE Ts2 offers brilliantly clear images, enabling more efficient cell culture observation.







ECLIPSE Ti2-U

Inverted Research Microscope with an excellent manual model

ECLIPSE Ti2-U provides an excellent base platform for accommodating a variety of research applications







The digital sight series and NIS-Elements are not for clinical diagnostic use

Specifications and equipment are subject to change without any notice or obligation on the part of the manufacturer. April 2021 ©2016-2021 NIKON CORPORATION

N.B. Export of the products* in this brochure is controlled under the Japanese Foreign Exchange and Foreign Trade Law. Appropriate export procedures shall be required in case of export from Japan.

*Products: Hardware and its technical information (including software)

⚠ WARNING

TO ENSURE CORRECT USAGE, READ THE CORRESPONDING MANUALS CAREFULLY BEFORE USING THE EQUIPMENT.



NIKON CORPORATION

Shinagawa Intercity Tower C, 2-15-3, Konan, Minato-ku, Tokyo 108-6290, Japan phone: +81-3-6433-3705 fax: +81-3-6433-3785

https://www.healthcare.nikon.com/en/

ISO 14001 Certified for NIKON CORPORATION

Nikon Instruments Inc. 1300 Walt Whitman Road, Melville, N.Y. 11747-3064, U.S.A.

phone: +1-631-547-8500; +1-800-52-NIKON (within the U.S.A. only) fax: +1-631-547-0299

https://www.microscope.healthcare.nikon.com/

Nikon Europe B.V. (headoffice)

Tripolis 100, Burgerweeshuispad 101, 1076 ER Amsterdam, Netherlands phone: +31-20-7099-000 fax: +31-20-7099-298

https://www.microscope.healthcare.nikon.com/en EU/

Nikon Instruments (Shanghai) Co., Ltd.

CHINA phone: +86-21-6841-2050 fax: +86-21-6841-2060 (Beijing branch) phone: +86-10-5831-2028 fax: +86-10-5831-2026 (Guangzhou branch) phone: +86-2-3882-0551 fax: +86-2-3882-0580 https://www.microscope.healthcare.nikon.com/zh_CN/

Nikon Canada Inc.

CANADA phone: +1-905-625-9910 fax: +1-905-602-9953

Nikon France, Succursale de Nikon Europe B.V. FRANCE phone: +33-1-4516-4516 fax: +33-1-4516-4505

Nikon Deutschland, Zweigniederlassung der Nikon Europe B.V.

GERMANY phone: +49-211-9414-888 fax: +49-211-9414-322

Nikon Italy, Branch of Nikon Europe B.V. ITALY phone: +39-055-300-9601 fax: +39-055-300-993

Nikon Europe B.V., Amsterdam, Zweigniederlassung Schweiz (Egg/ZH)

SWITZERLAND phone: +41-43-277-2867 fax: +41-43-277-2861

Nikon UK, Branch of Nikon Europe B.V.

UNITED KINGDOM phone: +44-208-247-1717 fax: +44-208-541-4584

Nikon Österreich, Zweigniederlassung der Nikon Europe B.V.

AUSTRIA phone: +43-1-972-6111 fax: +43-1-972-6111-40

Nikon Singapore Pte Ltd

SINGAPORE phone: +65-6559-3651 fax: +65-6559-3668 Nikon Instruments Korea Co., Ltd.



^{*2} NAMC (Nikon Advanced Modulation Contrast) is Nikon's unique modulation contrast observation method which provides stereoscopic images similar to DIC observation, even with samples on plastic dishes.

^{*3} Emboss contrast is Nikon's unique contrast observation method. It provides pseudo-three-dimensional images using focal illumination, which gives high contrast to samples.